

<Purification of VHH antibody by Protein A>

We investigate the ability of our Protein A resin product (Code; AAR-025) for purification of VHH antibody.

[Purification of VHH]

VHH antibody (+BSA) is applied to Protein A resin products (Our product, Competitor G and K). (*A; applied sample*)


↓ Unbound material (*F; flow-through fraction*)

wash by PBS

1

100mM Glysine, pH2.5 (E; elute fraction)

SDS-PAGE analysis (CBB stain)

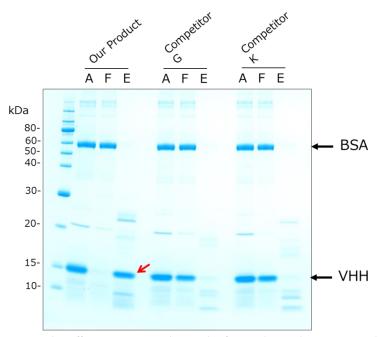

VHH antibody can be specifically captured and eluted in the use of our product for affinity resin, and not captured by competitor G, K products.

Reference)

Analytical Biochemistry 477 (2015) 92-94

Journal of Biotechnology 78 (2000) 11-21

Sales)

Please contact to Cosmo Bio (export@cosmobio.co.jp), our wholesaler.

